




High Power Crimping Tools

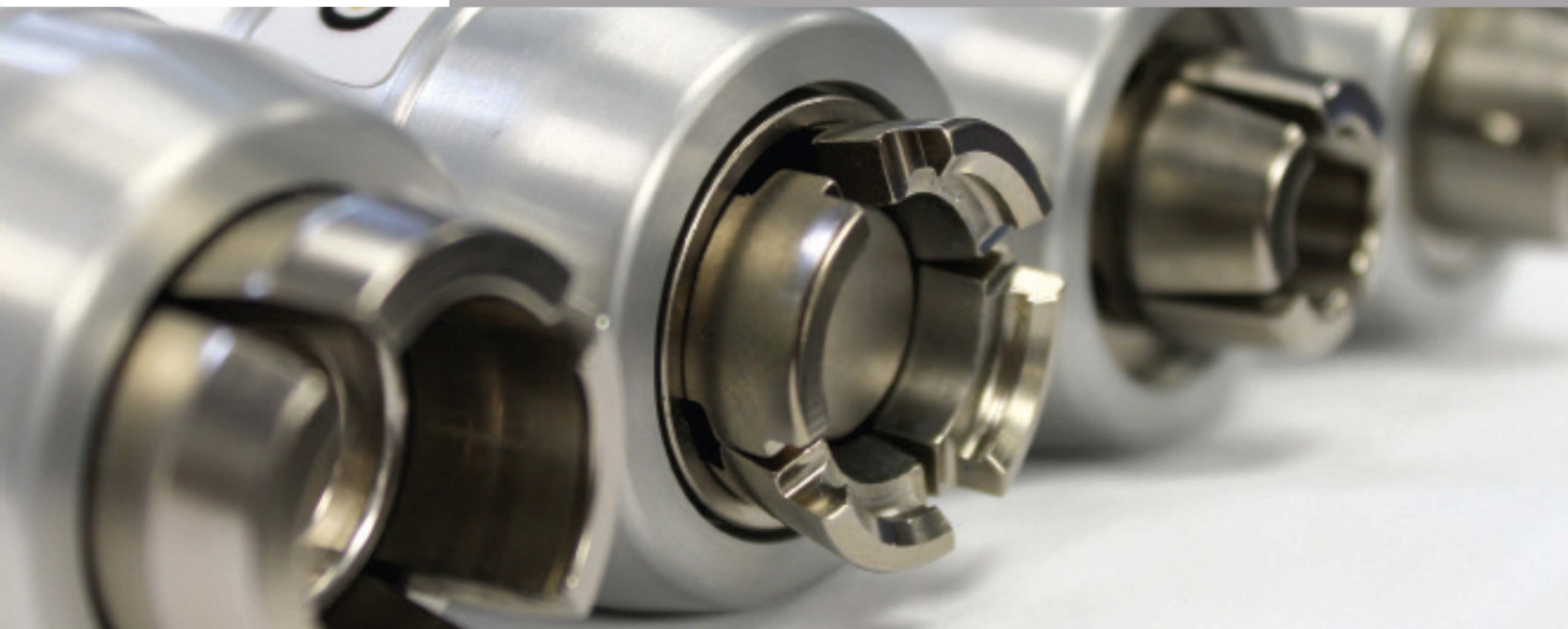

- High power version of electronic crimping tool has a fixed power supply and cord. (No battery)
- Uses interchangeable jaw sets
- OLED screen displays tool size, crimp setting and cycle result
- Settings menu shows lifetime statistics, recent cycle log, 8 language options and reset functions
- Strong enough for all steel and magnetic caps
- Convenient controls on top of unit to adjust crimp
- For flip-off caps, available in 13 mm and 20 mm sizes*
Not for use with Flip-up caps
- Optional base and mounting kit

The CS High Power Crimping Tool is designed with interchangeable jaws, so that you only need one tool for various functions in your laboratory. The jaw sets are available as crimper and decappers which fit standard lab vial seals with aluminum or two-part aluminum/steel caps. The tools are ergonomically designed to reduce strain and arm injury associated with many manual crimping tools. Jaw sets are available in 8 mm (crimper only), 11 mm, 13 mm, and 20 mm cap sizes, including flip-off caps in 13 mm and 20 mm sizes.

Switching Jaw Sets: Jaw sets are easy to interchange. Simply insert the jaw set into the bushing at the bottom of the tool. Push up and twist until the set locks into position. To remove, push the outside button and rotate.

Description	Item No.
High Power Electronic Crimping Tool	029150-V6
Base with Mounting Kit	029158
8 mm Crimper Jaw Set	029151
11 mm Crimper Jaw Set	029152
13 mm Crimper Jaw Set	029154
20 mm Crimper Jaw Set	029156

Description	Item No.
13 mm Jaw Set, for Flip Off Caps	029154-F0
20 mm Jaw Set, for Flip Off Caps	029156-F0
11 mm Decapper Jaw Set	029153
13 mm Decapper Jaw Set	029155
20 mm Decapper Jaw Set	029157